Product information

Information about other products is available at: www.demeditec.com

User's Manual

Pistachio ELISA

Enzyme Immunoassay for the quantitative determination of Pistachio in food

DEPISE01

96

Sensitivity (Pistachio)	0.13 ppm
Recovery	81-96 %
Incubation Time	60 min

1. GENERAL INFORMATION

Pistachio (*Pistacia vera*) belongs to the family of Anacardiaceae. With about 21 % the fraction of proteins in pistachio seed is very high. Some of these proteins, like the 2S albumin Pis v 1, the 11S globulin Pis v 2 or the 7S vicillin Pis v 3 are known for being allergenic. Many of them are heat resistant making them stable to different production processes. For this reason pistachio represents an important food allergen.

For pistachio-allergic persons hidden pistachio allergens in food are a critical problem. Already very low amounts of pistachio can cause allergic reactions, which may lead to anaphylactic shock in severe cases. Because of this, pistachio-allergic persons must strictly avoid the consumption of pistachio containing food. Cross-contamination, mostly in consequence of the production process, is often noticed. This explains why in many cases the existence of pistachio residues in food cannot be excluded. For this reason sensitive detection systems for pistachio residues in foodstuffs are required.

The **Pistachio ELISA** represents a highly sensitive detection system for pistachio and is particularly capable of the quantification of residues in cookies, cereals, ice-cream, chocolate and sausage.

2. PRINCIPLE OF THE TEST

The **Pistachio** quantitative test is based on the principle of the enzyme linked immunosorbent assay. An antibody directed against pistachio proteins is bound on the surface of a microtiter plate. Pistachio containing samples or standards are given into the wells of the microtiter plate. After 20 minutes incubation at room temperature, the wells are washed with diluted washing solution to remove unbound material. A peroxidase conjugated second antibody directed against pistachio proteins is given into the wells and after 20 minutes of incubation the plate is washed again. A substrate solution is added and incubated for 20 minutes, resulting in the development of a blue colour. The colour development is terminated by the addition of a stop solution, and the colour turns yellow. The yellow colour is measured photometrically at 450 nm. The concentration of pistachio is directly proportional to the colour intensity of the test sample.

3. PRECAUTIONS

Full compliance of the following good laboratory practices (GLP) will determine the reliability of the results:

1. Prior to beginning the assay procedure, bring all reagents to room temperature (20-25°C).
2. All reagents should be mixed by gentle inversion or swirling prior to use. Do not induce foaming.
3. Once the assay has been started, all subsequent steps should be completed without interruption and within the recommended time limits.
4. Replace caps in all the reagents immediately after use. Do not interchange vial stoppers.
5. Use a separate disposable tip for each specimen to prevent cross-contamination.
6. All specimens and standards should be run at the same time, so that all conditions of testing are the same.
7. Do not mix components from different batches.
8. Do not use reagents after expiration date.
9. Check both precision and accuracy of the laboratory equipment used during the procedure (micro-pipets, ELISA reader etc.).

4. HEALTH AND SAFETY INSTRUCTIONS

1. Do not smoke or eat or drink or pipet by mouth in the laboratory.
2. Wear disposable gloves whenever handling patient specimens.
3. Avoid contact of substrate and stop solution with skin and mucosa (possible irritation, burn or toxicity hazard). In case of contact, rinse the affected zone with plenty of water.
4. Handling and disposal of chemical products must be done according to good laboratory practices (GLP).

5. REAGENTS

The kit contains reagents for 96 determinations. They have to be stored at 2-8°C. Expiry data are printed on the labels of the bottles and the outer package.

1. Microtiter plate consisting of 12 strips with 8 breakable wells each, coated with anti-pistachio antibodies.
2. Pistachio Standards (0; 1; 4; 10; 40 ppm of pistachio): 5 vials with 1.0 mL each, dyed red, ready-to-use.
3. Conjugate (anti-pistachio-peroxidase): 15 mL, dyed red, ready-to-use.
4. Substrate Solution (TMB): 15 mL, ready-to-use.
5. Stop Solution (1 N acidic solution): 15 mL, ready-to-use.
6. Extraction and sample dilution buffer (Tris): 2 x 120 mL as 10x concentrate, dyed red. Dilute 1+9 with distilled water. Stored at 4°C the diluted buffer is stable for at least one week. If during the cold storage crystals precipitate, the concentrate should be warmed up to 37°C for 15 minutes.
7. Washing Solution (PBS + Tween 20): 60 mL as 10x concentrate. Dilute 1+9 with distilled water. Stored at 4°C the diluted buffer is stable for at least 4 weeks. If during the cold storage crystals precipitate, the concentrate should be warmed up to 37°C for 15 minutes.
8. Plastic bag to store unused microtiter strips.
9. Instruction Manual.

6. ADDITIONAL INSTRUMENTATION AND REAGENTS (NOT PROVIDED)

Instrumentation

- 100 - 1000 µL micropipets
- Volumetric flask
- Analytical balance
- Mortar, mixer
- Water bath
- Centrifuge
- ELISA reader (450 nm)

Reagents

- double distilled water

7. SAMPLE PREPARATION

Due to high risk of cross-contamination all applied instruments like applicator, mortar, glass vials etc. have to be **cleaned thoroughly** before and after each sample. Pistachio proteins could adhere to different surfaces. To identify possible cross-contamination caused by previous extractions it is strongly recommended to note the sequence of the extractions.

The following sample preparation should be applied for solid samples:

1. To maximize homogeneity and representativeness of the sample drawing, a minimum of 5 g sample should be pulverized finely in a mortar, impact mill etc.
2. 1 g of the homogenized mixture is suspended in 20 mL of **pre-diluted** extraction and sample dilution buffer. Afterwards the suspension is incubated for 15 min in a preheated water bath at 60°C. To ensure good homogeneity, the samples should be shaken every two minutes.
3. The samples are centrifuged for 10 minutes at 2000 g. If it is not possible to separate the supernatant from the precipitate completely, the suspension should be filtrated if necessary.
4. 100 µL of particle-free solution are applied per well. If the results of a sample are out of the measuring range, further dilution with the **pre-diluted** extraction and sample dilution buffer is necessary. The additional dilution has to be considered when calculating the concentration.

The following sample preparation should be applied for liquid samples:

1 mL of liquid sample is diluted in 19 mL of **pre-diluted** extraction and sample dilution buffer. Afterwards the suspension is incubated for 15 min in a preheated water bath at 60°C. To ensure good homogeneity, the samples should be shaken every two minutes. The process is continued at point 3 of solid sample extraction process.

8. PROCEDURE

The washing solution is supplied as 10x concentrate and has to be **diluted** 1+9 with double distilled water before use.

In any case the **ready-to-use** standards provided should be determined twofold. When samples in great quantities are determined, the standards should be pipetted once before the samples and once after the samples. For final interpretation the arithmetic mean is used for calculation.

In consideration of GLP and quality control requirements a duplicate measurement of samples is recommended.

The procedure is according to the following scheme:

1. Prepare samples as described above.
2. Pipet 100 µL **ready-to-use** standards or prepared samples in duplicate into the appropriate wells of the microtiter plate.
3. Incubate for 20 minutes at room temperature.
4. Wash the plate three times as follows: Discard the contents of the wells (dump or aspirate). Pipet 300 µL of diluted washing solution into each well. After the third repetition empty the wells again and remove residual liquid by striking the plate against a paper towel. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbencies.
5. Pipet 100 µL of conjugate (anti-pistachio-peroxidase) into each well.
6. Incubate for 20 minutes at room temperature.
7. Wash the plate as outlined in 4.
8. Pipet 100 µL of substrate solution into each well.
9. Allow the reaction to develop in the dark (e.g. cupboard or drawer; the chromogen is light-sensitive) for 20 minutes at room temperature.
10. Stop enzyme reaction by adding 100 µL of stop solution (1 N acidic solution) into each well. The blue colour will turn yellow upon addition.
11. After thorough mixing, measure absorbance at 450 nm (reference wavelength 620 nm), using an ELISA reader. The colour is stable for 30 minutes.

9. CALCULATION OF RESULTS

The ready-to-use standards are prepared for a direct determination of sample concentrations. The dilution of samples in the extraction process as described in the above stated sample preparation procedure is already considered. Additional dilution due to high sample concentration has to be accounted for.

1. Calculate the average optical density (OD 450 nm) for each set of reference standards or samples.
2. Construct a standard curve by plotting the mean optical density obtained for each reference standard against its concentration in ppm on semi-log graph paper with the optical density on the vertical (y) axis and the concentration on the horizontal (x) axis. Alternatively the evaluation can be carried out by software. In this case the 4-parameter method should be preferred.
3. Using the mean optical density value for each sample, determine the corresponding concentration of pistachio in ppm from the standard curve. Depending on experience and/or the availability of computer capability, other methods of data reduction may be employed.

10. TYPICAL STANDARD VALUES

The following table contains an example for a typical standard curve. The binding is calculated as percent of the absorption of the 40 ppm standard. These values are only an example and should not be used instead of the standard curve which has to be measured in each new test.

Pistachio (ppm)	% binding of 40 ppm
40	100
10	54
4	29
1	10
0	3

11. PERFORMANCE

Sensitivity

The limit of detection (LOD) of the **Pistachio test** is 0.13 ppm for the standard curve. Validation experiments with common matrices resulted in the following LODs [ppm].

Cookies	0.18
Cereals	0.01
Ice-cream	0.20
Chocolate	0.09
Sausage	0.17

The limit of quantification (LOQ) of the **Pistachio test** is 1 ppm.

Due to the variety of sample matrices and their influence on the blank, results less than the LOQ should be treated as negative.

Precision

Intra-assay Precision	6-8 %
Inter-assay Precision	5-12 %

Linearity

The serial dilution of spiked samples (cookies, cereals, ice-cream, chocolate and sausage) resulted in a dilution linearity of 90–116 %.

Cross-reactivity

For the following foods no cross-reactivity could be detected:

Almond	Coconut	Plum
Apricot	Cod	Poppy seed
Barley	Corn	Pork meat
Bean	Cow's milk	Potato
Beef	Egg	Pumpkin seed
Bovine gelatin	Guar gum	Rice
Brazil nut	Kiwi	Rye
Carob gum	Lentil	Saccharose
Carrot	Lupin	Sesame
Celery	Macadamia	Shrimp
Cherry	Mustard	Soy
Chestnut	Oats	Soy lecithin
Chicken meat	Pea	Tomato
Chickpea	Peanut	Wheat
Cocoa	Pine seed	

The following cross reactions were determined:

Cashew	12 %
Hazelnut	0.17 %
Pecan	0.0005 %
Sunflower seed	0.0002 %
Walnut	0.0008 %

Recovery

Mean recovery was determined by spiking samples with different amounts of pistachio:

Cookies	94 %
Cereals	96 %
Ice-cream	89 %
Chocolate	81 %
Sausage	90 %

12. REFERENCES

1. Noorbakhsh R, et al. (2011) – Pistachio allergy-prevalence and in-vitro cross-reactivity with other nuts. *Allergol Int*, 60(4):425-432
2. Lim PW (2010) – Development of an enzyme-linked immunosorbent assay for the detection of pistachio residues in processed foods. Dissertation, University of Nebraska-Lincoln
3. Noorbakhsh R, et al. (2010) – Influence of processing on the allergenic properties of pistachio nut assessed in vitro. *J Agric Food Chem*, 58:10231-10235
4. Ahn K, et al. (2009) – Identification of two pistachio allergens, Pis v 1 and Pis v 2, belonging to the 2S albumin and 11S globulin family. *Clin Exp Allergy*, 39(6):926-934

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